

Instruction Manual

Product	Size	Catalog Number
Cryo-SFM	30 ml	C-29910
	125 ml	C-29912

Product Description

PromoCell Cryo-SFM is a protein-free, chemically defined and animal component-free cryopreservation medium. The optimized formulation is based on methylcellulose, DMSO and other cryoprotectants. Cryo-SFM provides excellent results with all types of cells including primary human cells, stem cells and cell lines. Cells frozen in Cryo-SFM exhibit superior viability, attachment and subsequent growth performance after thawing.

Instructions for Use

Detach adherent cells using the PromoCell DetachKit (C-41200, C-41210, C-41220).

Resuspend the cells in cold or room temperature Cryo-SFM at a concentration of 0.1 – 10 million cells/ml. Transfer the cell suspension into cryo-vials, freeze them gradually (-1°C/minute) to -70°C or lower overnight and subsequently store the vials in liquid nitrogen.

Note: Each freeze/thaw cycle leads to a loss in proliferation potential and cell number, particularly when working with fragile primary cells. If freezing the cells, please do so at as low a passage number as possible.

Storage and Stability

Store Cryo-SFM at 4 to 8°C in the dark immediately after arrival. If stored properly, the product is stable until the

expiry date stated on the label.

Quality Control

All lots of PromoCell Cryo-SFM are subjected to comprehensive quality control tests. Each lot is routinely tested for its performance with primary human cells, for the absence of cytotoxicity, and for physical parameters such as osmolality and pH level. Approved in-house lots are used as a reference.

In addition, all lots of Cryo-SFM have been tested for the absence of microbial contaminants (fungi, bacteria).

Intended Use

The products are for *in vitro* research use only and not for diagnostic or therapeutic procedures. For safety precautions please see appropriate MSDS.

PromoCell GmbH

Sickingenstr. 63/65
69126 Heidelberg
Germany

Email: info@promocell.com
www.promocell.com

USA/Canada

Phone: 1 – 866 – 251 – 2860 (toll free)
Fax: 1 – 866 – 827 – 9219 (toll free)

Deutschland

Telefon: 0800 – 776 66 23 (gebührenfrei)
Fax: 0800 – 100 83 06 (gebührenfrei)

France

Téléphone: 0800 90 93 32 (ligne verte)
Téléfax: 0800 90 27 36 (ligne verte)

United Kingdom

Phone: 0800 – 96 03 33 (toll free)
Fax: 0800 – 169 85 54 (toll free)

Other Countries

Phone: +49 6221 – 649 34 0
Fax: +49 6221 – 649 34 40

© PromoCell GmbH